

# SILMARK CART

*Lifting fingermarks from irregular or textured surfaces  
Available in Black, White, and Clear*

FINGERMARK ON POLYSTYRENE SURFACE, DEVELOPED WITH MAGNETIC JET BLACK FINGERPRINT POWDER. LIFTED WITH SILMARK CART WHITE WITH ATTACHED SILICONE PAPER (CAT.NO. C-106010).



Silmark CART is a 2-part addition curing silicone compound in a cartridge. Silmark CART is available in various colours: gray and brown for toolmark casting and black, white, and clear for lifting of powdered fingermarks.

Using the dispensing gun the compound is easily applied to textured surfaces like Hammerite painted metal and polystyrene packing materials.

There is no need to manually mix the two components, this is done by the static mixing tip which is mounted on the cartridge prior to each use. With the optional spreader tip attached to the static mixer a 25 mm/1 inch wide strip of silicone material is applied over the fingermark to be lifted.

A sheet of silicone paper can be applied to the silicone material shortly after applying it to the surface containing the fingermark. Gently rubbing the back of the paper ensures the silicone material is fully pressed into the texture of the surface.

This will not only assist in getting the powdered fingermark off the surface as complete as possible, in many cases it will also provide a flat backside of the lift which makes it easier to handle during photography. The paper adheres to the silicone lift and offers ample space to make notes on.

Setting times vary with temperature, with ambient and material temperatures of 20°C/68°C this is around 40 seconds and some 80 seconds at 10°C/50°F.



**BVDA**  
FORENSIC PRODUCTS

BVDA INTERNATIONAL BV • INFO@BVDA.NL • WWW.BVDA.COM



DISPENSING GUN FOR 50 ML CARTRIDGES (CAT.NO. C-8210).



WITH A SPREADER TIP (CAT.NO. C-8350) ATTACHED TO THE STATIC MIXER A STRIP OF 25 MM WIDE MATERIAL IS SIMPLY APPLIED TO THE FINGERMARK.



LIFT OF A FINGERMARK DEVELOPED WITH BLACK FINGERPRINT POWDER USING SILMARK CART CLEAR (CAT.NO. C-2505).



## SILMARK CART FOR FINGERMARK LIFTING

CAT.NO.	DESCRIPTION
C-1405	Silmark CART White, 75 ml
C-2105	Silmark CART Black, 75 ml
C-2505	Silmark CART Clear, 75 ml

## AUXILIARY PRODUCTS

CAT.NO.	DESCRIPTION
C-8200	Silmark CART dispensing gun for 50 and 75 ml cartridges
C-8210	Silmark CART dispensing gun for 50 ml cartridges only
C-8300	Static mixing tips, 50/p
C-8350	Spreader tips 25 mm/1", 50/p
C-106010	Siliconized paper (one side only) 10 x 14 cm / 4 x 5,6", 100/p

## SILMARK CART FOR TOOLMARK CASTING

CAT.NO.	DESCRIPTION
C-1105	Silmark CART Grey, 75 ml
C-1106	Silmark CART Grey, 50 ml
C-1005	Silmark CART Wintry Grey, 75 ml
C-1006	Silmark CART Wintry Grey, 50 ml
C-2405	Silmark CART Brown, 75 ml
C-2406	Silmark CART Brown, 50 ml
C-2305	Silmark CART Wintry Brown, 75 ml
C-2306	Silmark CART Wintry Brown, 50 ml